

# African American Admixture Panel

Targeted multi-sample genotyping panel from Illumina delivers high-quality data and is a powerful tool for conducting admixture mapping studies.

## INTRODUCTION

Admixture mapping can be a powerful method for identifying genomic regions harboring common disease susceptibility loci in recently mixed populations. The Illumina African American Admixture Panel contains 1,509 SNPs for admixture mapping studies in African American populations. This specialized content was developed and validated in collaboration with Dr. David Reich at Harvard Medical School and the Broad Institute of MIT and Harvard.

Admixture mapping requires a genome-wide panel of markers that can differentiate chromosomal segments in admixed individuals by their ancestral origins. African American populations are of particular interest for admixture mapping studies due to the recent mixture of African and European populations. There have been an average of six generations since African and European populations intermixed after relocation into North America. Thus, there has been relatively little

### HIGHLIGHTS OF THE AFRICAN AMERICAN ADMIXTURE PANEL

- **High-Quality Data:** Proven GoldenGate® technology
- **Targeted Content:** Powerful method to detect disease susceptibility loci
- **High Throughput:** Quick and efficient assay of 96 samples simultaneously

recombination between African- and European-derived chromosomes in African American populations. This low amount of recombination has yielded chromosomal segments of either European or African ancestry that are typically tens of megabases long (average ~17cM<sup>1</sup>). The Illumina African American Admixture Panel enables a highly-efficient scan across the entire genome for genetic disease risk factors using a minimum number of markers. By targeting markers to these large single-ancestry derived chromosomal segments, only relevant variance in the admixed population is interrogated. By specifically targeting variation between European and African chromosomes, relatively few markers provide the same power as hundreds-fold more randomly

chosen SNPs. Consequently, many more samples can be analyzed with an intelligently targeted genotyping panel at much lower cost.

### ADMIXTURE MAPPING WITH A TARGETED GENOTYPING PANEL

The Illumina African American Admixture Panel facilitates detection of susceptibility loci of genetic variants that differ in frequency between ancestral African and European populations. Population admixture events between previously geographically separated populations, such as occurred between African and European populations in America, introduce new linkage disequilibrium (LD). Samples from disease cases can be genotyped and compared to find regions of the genome where the proportion of

FIGURE 1: AFRICAN AMERICAN ADMIXTURE PANELS ARE DEPLOYED IN MULTI-SAMPLE SENTRIX ARRAY MATRIX FORMAT




Admixture mapping experiments are facilitated by high-throughput sample analysis. Marker panels are deployed on a 96-sample Sentry® Array Matrix.

African ancestry is different from the genome-wide average. Regions with African ancestry deviating from average are likely to contain genetic variants associated with disease. Finding a disease-associated region with this panel allows more in-depth focused studies aimed at uncovering direct causes. Illumina iSelect™ Infinium® Custom Genotyping Panels and sequencing using Illumina Genome Analyzer powered by Solexa® Technology are both excellent options for subsequent fine-mapping.

Association studies using Illumina panels based on GoldenGate Assay technology allow for broad genomic screening, and are highly efficient due to the rationally targeted SNP marker selection. A streamlined panel of 1,509 targeted loci yields approximately 75% of the power for admixture mapping, as would be expected with 300,000–1,000,000 loci required for dense non-specific whole-genome scans. Furthermore, efficiency is greatly enhanced by using the high-throughput multi-sample Sentrix Array Matrix format, which permits 96 simultaneous assays.

Admixture mapping is a useful method to study diseases that differ in frequency between different populations. Diseases where African American admixture mapping is anticipated to be effective include

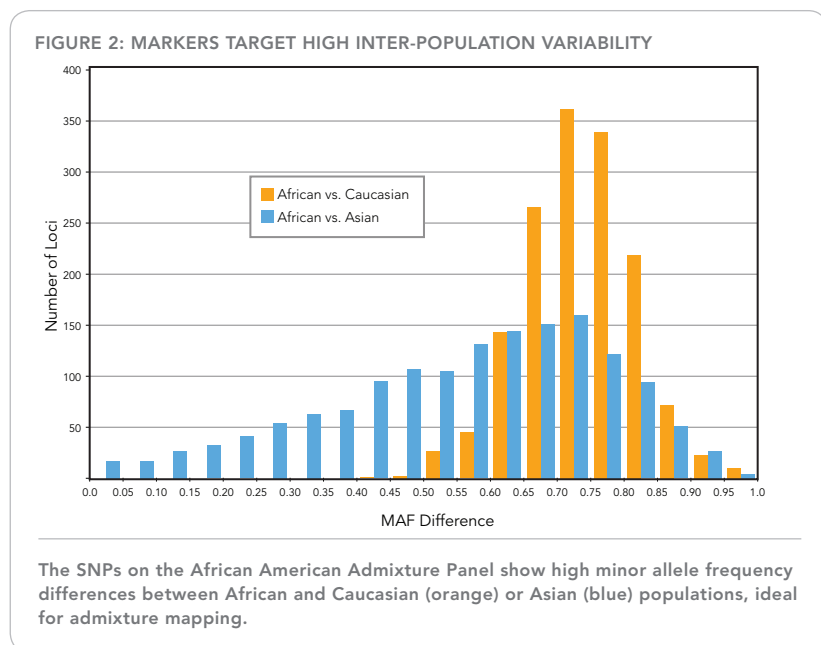

prostate cancer<sup>2,3</sup>, Type-II diabetes, autoimmune diseases (e.g. multiple sclerosis<sup>4</sup>), hypertension<sup>5</sup>, asthma<sup>6</sup>, and obesity<sup>7</sup>. Earlier versions of the panel have allowed researchers to successfully localize genes associated with multiple sclerosis<sup>4</sup>, prostate cancer<sup>2</sup>, and serum markers of inflammation<sup>8</sup>. For example, a strong association between a 3.8Mb chromosomal region (8q24) and susceptibility to prostate cancer in African Americans was recently found using an earlier version of this panel<sup>2</sup>.

#### TARGETED CONTENT

The optimal set of loci for admixture mapping are those with high frequency differences between African and European populations (Figure 2). This African American Admixture Panel was designed using a multi-step selection process. Millions of SNP loci across the genome were screened bioinformatically to determine those most likely to provide ancestry information in African American populations<sup>8</sup>. The resulting thousands of SNP loci in several African and European populations were then screened and verified using the Illumina GoldenGate Assay. SNPs

**TABLE 1: AFRICAN AMERICAN ADMIXTURE PANEL DATA QUALITY (285 SAMPLES, 15 REPLICATES, 60 TRIPLICATES)**

POPULATION	AVG MAF (LOCI)	CALL RATE	AVG HET (DNA)	REPRODUCIBILITY	HERITABILITY
Caucasian (CEU)	11.92%	99.89%	19.46%		
Asian (JPT+CHB)	23.49%	99.92%	30.66%		
African (YOR)	14.32%	99.86%	22.87%		
Overall		99.89%		100%	100%

failing Hardy-Weinberg equilibrium testing ( $P > 0.01$ ) were eliminated. This optimized set of markers detects a majority of the variation differentiating African and European chromosomal segments.

#### HIGH-QUALITY DATA

The African American Admixture Panel was subjected to rigorous functional testing to ensure high-quality, reproducible data. Samples derived from each of three HapMap population groups (Caucasian, Asian, and

African) exhibited call rates greater than 99.8%, with 100% reproducibility across 15 replicates and 60 triplicates (Table 1).

#### ILLUMINA SOLUTIONS FOR GENOTYPING

The high quality and low cost per genotype of the African American Admixture Panel are made possible by powerful Illumina technologies that include the GoldenGate Assay with multi-sample Sentrix Array Matrix format. Illumina genotyping

solutions enable linkage and association mapping, as well as high-resolution, genome-wide scans of greater than 1,000,000 markers per BeadChip. Illumina genotyping panels with standard or custom content can be accessed via FastTrack Genotyping Services or with an Illumina System. The combination of Illumina's proprietary assay technologies and intelligent content selection delivers the most comprehensive DNA analysis solutions now available.

#### ORDERING INFORMATION

CATALOG NO.	PRODUCT	DESCRIPTION
GT-17-230	African American Admixture Panel	One oligo pool (OPA) for 1,509 SNP loci. Sufficient for 96 DNA samples.
GT-95-201	Single Use Activation Kit (576 Samples)	Used in combination with the GoldenGate Assay Kit. Contains reagents for six 96-well plates.
GT-95-205	GoldenGate Assay Kit II with UDG (96 Samples)	Sufficient reagent for preparing genotyping reactions for 96 DNA samples. Contains UDG enzymes for contamination control.
FA-12-112	Sentrix Universal-96 Array Matrix (1536-plex)	A single Sentrix Array Matrix that can process 96 samples with up to 1536 assays/sample.

## REFERENCES

- (1) Patterson N, Hattangadi N, Lane B, Lohmueller KE, Hafler DA et al. (2004) Methods for high-density admixture mapping of disease genes. *Am J Hum Genet* 74: 979-1000.
- (2) Freedman ML, Haiman CA, Patterson N, McDonald GJ, Tandon A et al. (2006) Admixture mapping identifies 8q24 as a prostate cancer risk locus in African-American men. *Proc Natl Acad Sci U S A* 103: 14068-14073.
- (3) Haiman CA, Patterson N, Freedman ML, Myers SR, Pike MC et al. (2007) Multiple regions within 8q24 independently affect risk for prostate cancer. *Nat Genet* 39: 638-644.
- (4) Reich D, Patterson N, De Jager PL, McDonald GJ, Waliszewska A et al. (2005) A whole-genome admixture scan finds a candidate locus for multiple sclerosis susceptibility. *Nat Genet* 37: 1113-1118.
- (5) Zhu X, Luke A, Cooper RS, Quertermous T, Hanis C et al. (2005) Admixture mapping for hypertension loci with genome-scan markers. *Nat Genet* 37: 177-181.
- (6) Barnes KC, Grant AV, Hansel NN, Gao P, Dunston GM (2007) African Americans with asthma: genetic insights. *Proc Am Thorac Soc* 4: 58-68.
- (7) Cossrow N, Falkner B (2004) Race/ethnic issues in obesity and obesity-related comorbidities. *J Clin Endocrinol Metab* 89: 2590-2594.
- (8) Smith MW, Patterson N, Lautenberger JA, Truelove AL, McDonald GJ et al. (2004) A high-density admixture map for disease gene discovery in African Americans. *Am J Hum Genet* 74: 1001-1013.
- (9) Reich D, Patterson N, Ramesh V, De Jager PL, McDonald GJ et al. (2007) Admixture mapping of an allele affecting interleukin 6 soluble receptor and interleukin 6 levels. *Am J Hum Genet* 80: 716-726.

## ADDITIONAL INFORMATION

Visit our website or contact us at the address below to learn more about Illumina DNA Analysis solutions.

## Illumina, Inc.

## Customer Solutions

9885 Towne Centre Drive  
 San Diego, CA 92121-1975  
 1.800.809.4566 (toll free)  
 1.858.202.4566 (outside the U.S.)  
[techsupport@illumina.com](mailto:techsupport@illumina.com)  
[www.illumina.com](http://www.illumina.com)

## FOR RESEARCH USE ONLY

© 2007 Illumina, Inc. All rights reserved.

Illumina, Solexa, Making Sense Out of Life, Oligator, Sentrix, GoldenGate, DASL, BeadArray, Array of Arrays, Infinium, BeadXpress, VeraCode, IntelliHyb, iSelect, and CSPro are registered trademarks or trademarks of Illumina. All other brands and names contained herein are the property of their respective owners.  
 Pub. No. 370-2007-011 12Jul07

